



530 Main Street, Acton, MA 01720

**Phone:** (978)263-3584, **Fax:** (978)263-5086

**Web Site:** [www.acton-research.com](http://www.acton-research.com)

## **Supplemental Instructions for External Shutter**

Per Customer specifications this SpectraPro is supplied with an External Shutter to mount on the outside of any standard SpectraPro entrance slit assembly. To attach the External Shutter to the slit assembly, insert the supplied #8-32 socket head cap screws into the counter bored holes in the Shutter and screw the Shutter fast to the slit, assuring that the label “TOP” is on the top. The shutter is normally closed and needs to be activated before light will pass through the instrument.